**IKA®** RH basic 2  
**IKA®** RH basic KT/C  
**IKA®** RH digital KT/C

IKA RH basic 2

IKA RH basic KT/C

IKA RH digital KT/C

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## CE-DECLARATION OF CONFORMITY

**GB**

We declare under our sole responsibility that this product corresponds to the regulations 89/336/EEC; 92/31/EEC and 73/23EEC and conforms with the standards or standardized documents EN 61010-1 (IEC/CEI 61010-1) and EN 61326-1 (IEC 61326-1).

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## Warranty

You have purchased an original IKA laboratory machine which meets the highest engineering and quality standards.

In accordance with IKA guarantee conditions, the guarantee period is 24 months. For claims under the guarantee please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

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## Contens

	Page
Warranty	2
Safety instructions	11
Proper use	12
Unpacking	12
Commisioning	12
Starting	12
Function stirring	13
Function heating	13
Limiter of safety temperature	14
Adjusting the Safety Temperature limit "Safe Temp" (Vers. KT)	14
Controlling the Medium Temperature via ETC 1 or Contact Thermometer (Vers. KT)	15
Error messages	16
Mounting - Support H 16 V	16
Maintenance and cleaning	16
Accessories	17
Technical data	18
Listofspare parts RH basic 2 / RH basic KT/C / RH digital KT/C	28
Spare parts diagram RH basic 2 / RH basic KT/C / RH digital KT/C	29-31

## Safety instructions

The IKA RH basic 2; RH basic KT/C or RH digital KT/C was not constructed for operations in dangerous atmospheres, for mixing dangerous substances or for operation under water. The instrument may not be operated in rooms with explosion hazard. Furthermore, it has to be placed upon a fire-proof and/or not burnable surface.

**Attention - Magnetism!** Effects of the magnetic field have to be taken into account (e.g. data carriers, cardiac pacemakers ...). Please make sure that the mains cable does not contact the heating plate. The power cable must be replaced by a cable of equal quality and capacity.

**Version RH basic KT/C and RH digital KT/C:** Examine regularly (we recommend every 4 weeks) the function of the safety temperature limiter. For this the safety temperature must be changed around at least 1/4 revolution of the value already stopped and be accomplished the attitude as under "adjusting the safety temperature delimitation SAFE temp" described. By the exceeding of the adjusted safety temperature (e.g. rotary button "temp" posed on right end stop) you can examine the function. The "temp" pilot lamp must expire and the heating must be lasting switched off.

When using PTFE-coated magnetic bars, the following have to be noted: *Chemical reactions of PTFE occur in contact with molten or dissolved alkaline and alkaline - earth metals, as well as with fine-particled powders of metals of the 2. and 3. group of the periodical system at temperatures above 300-400°C. Only elementary fluorine, chlorine trifluoride und alkaline metals do attack PTFE, halogen hydrocarbons have a reversibly swelling effect.*

Source: Römpps Chemie-Lexikon and „Ullmann“ Bd.19

Exercise caution when touching the housing parts and the heating plate!

**Risk of burns!** The heating plate can reach temperatures in excess of 300°C.

Please consider possible endangering due to release of toxic or flammable gas caused by heated medium.

Before placing the device in service, please ensure that the turn dial for adjusting the speed is against the left stop, since the device will begin to run at the last speed that was set.

After an interruption in the power supply during an agitation procedure, the device will start running again by itself.

Because of increased temperature of the housing, you should only move the unit when it has been turned off and has cooled down.

The instrument may not be transported on a H 16 V support rod or on the connection line.

Clicking noises may occur as a result of material expansion during heating and cooling.

**Caution:** Only process and heat up any media that has a flash point higher than the adjusted safety temperature limit (max 400°C for RH basic 2, 100°C to max. 400°C for RH basic KT/C and RH digital KT/C).

Please observe the relevant safety information and guidelines as well as work protection and accident prevention requirements for use in the laboratory.

Only a specialist may open the device.

Use only original IKA accessories.

## Proper use

The IKA RH basic 2; RH KT; RH basic KT/C or the IKA RH digital KT/C is a stirring instrument with heating function. It is used in e.g. laboratories of the chemical industry, in schools as well as in pharmacies. The instrument is suitable for tempering substances which are placed on the heating plate in vessels. By means of a built-in stirring drive and with the help of a magnetic bar within the vessel, the substances can be stirred at the same time. The mixing intensity is dependent on the motor speed and on the size of the magnetic bar. The Der RH basic KT/C and RH digital KT/C also has a contact thermometer connection in compliance with requirements of DIN 12 878 and an adjustable safety circuit.

## Unpacking

Please unpack the equipment carefully and check for any damages. It is important that any damages which may have arisen

during transport are ascertained when unpacking. If applicable a fact report must be set immediately (post, rail or forwarder).

The delivery scope covers: an RH basic 2 / RH basic KT/C or RH digital KT/C and operating instructions.

## Commissioning

Check whether the voltage specified on the type plate matches the mains voltage available. The power socket used must be earthed (protective earth conductor contact). Before tuning on the machine, adjust the turn dial for controlling the speed so that it is against the left stop. If these conditions are met, the device is ready to operate after plugging in the mains plug. If these procedures are not followed, safe operation cannot be guaranteed and/or the equipment may be damaged.

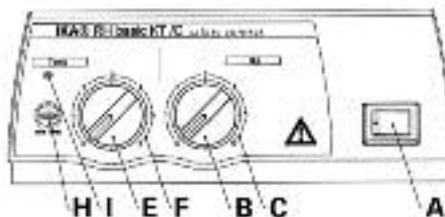
Observe the ambient conditions (temperature, humidity, etc.) listed under Technical Data.

## Starting


The heating stirring features can be turned on and off by pressing the unit switch **A**.

**Version RH basic 2 and RH basic KT/C:** The ready for operation and turned on states of the device are indicated by the

green lamp in the unit switch lighting up.

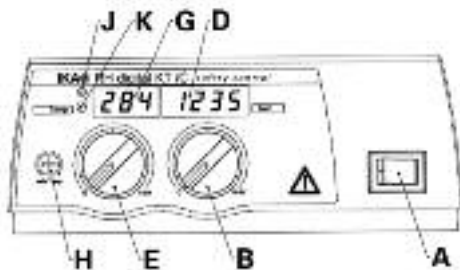**Version RH digital KT/C:** When you switch on the stirrer it will run a self-test and all the LED segments of the displays will light

up for a few seconds. After this, the LED's (**D**; **G**) light up to indicate that the unit is powered up and ready for operation.

## Function Stirring

The stirring rod in the glass beaker with a maximum length of 60mm is driven by a permanent magnet located directly under the placing and heating surface. It allows magnetic field lines to pass through it. The permanent magnet is fastened directly onto the drive shaft of the motor.

The actual speed depends on the load and voltage. Note that fluctuations in the mains power voltage within the allowable tolerance and changes in viscosity of the materials to be stirred resulting from processes they undergo may also cause minor fluctuations in the speed.

**Version RH basic 2 and RH basic KT/C:** The speed of the motor can be adjusted on the "Mot" dial (**B**) with a scale from 0 to 6 (**C**). The scale values from 0 to 6 on this dial correspond roughly to speeds ranging from 0 to 2000 rpm. The maximum speed of 2000 rpm is reached when the dial switch is turned all the way to the right and there is no load on the motor.

**Version RH digital KT/C:** Set the motor speed using the "Mot" knob (**B**); the knob has a graduated scale from to maximum rpm.

The motor speed is shown on the LED display (**D**). When the motor is stopped the speed "0" will be displayed. When the motor is operating, the display will show the current speed in the range 100 to 2000 rpm.

## Function Heating

The temperature of the heating plate of the instrument is kept constant by a control circuit. In addition, it is monitored by a safety circuit. The required temperature sensor (a thermo-couple) is installed in the heating plate. The **RH basic** unit has a stainless steel surface plate with a heating output of 400Watt.

**RH basic KT/C and RH digital KT/C** units have a enamel-coated stainless steel surface plate with a heating output of 500 Watt.

**Version RH basic und RH basic KT/C:** The temperature of the heating plate can be adjusted on the „Temp“ turn dial (**E**) with a corresponding scale from 0 to 6 (**F**). The scale values from 0 to 6 correspond approximately to area temperatures up to a max. of 320 °C. When the setting is at 0, the device does not heat the medium up at normal room temperature.

During the heating phase, the signal light (**I**) lights up orange (orange = power is being supplied to the heating plate; signal light off = no power being supplied)

**Version RH digital KT/C:** The temperature of the heating plate can be adjusted on the „Temp“ turn dial. When you change the knob setting, the LED display (**G**) will show the SETPOINT temperature and the warning light (**K**) will light up. After approx. 3 seconds the LED display will switch back to showing the ACTUAL temperature and the warning light (**K**) will switch off.

The temperature indication now constantly changes between actual temperature (25 sec) and target temperature (5 sec).

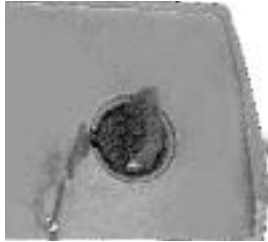
When you select the SETPOINT temperature and as the unit heats up, the warning light (**K**) will remain off and the LED display (**G**) will show the ACTUAL temperature of the surface plate in the range from 0°C to +320°C.

The red warning light (**J**) will light up while heating is in progress. (Red warning light ON = surface plate heating power supply is ON; red warning light OFF = surface plate heating power is OFF).

#### PLEASE NOTE!

**The heating plate of the version RH basic KT/C and RH digital KT/C only heats up, if the contact plug has been plugged** (the contact plug is undetachably fitted to the instrument), or if an **ETC 1** and/or a **contact thermometer** has been connected.

## Limiter of safety temperature

The safety temperature limiter makes it possible to adjust the maximum temperature of the heating plate (safety temperature). This ensures that the temperature of heating plate does not exceed the flame point of the material. The value to be set must be determined by trial and error and should generally fall about 20K below the flame point of the medium.

RH basic 2 devices have a safety circuit with a fixed setting that limits the temperature of a heating plate to 400°C.

RH basic KT/C and RH digital KT/C devices have an adjustable safety circuit. The maximum temperature of the heating plate can be adjusted within a range from 100°C to max. 400°C.

## Adjusting the safety temperature limit "Safe Temp" (Vers. KT)

**Version RH basic KT/C and RH digital KT/C:** RH basic KT/C and RH digital KT/C versions: The "Safe Temp" set screw (**H**) is to the left of the "Temp" knob (**E**) on the front operator panel.

You can use the „Safe Temp“ set screw (**H**) of the safety temperature limiter to adjust the desired safety temperature. If the desired safety temperature is exceeded by a higher adjusted target temperature, or if there is a defect in the instrument controller, the heating turns off permanently in this state and **the stirring function continues to run at the speed that was set prior to the malfunction.**

- Using a straight-edge screwdriver, turn the "Safe Temp" adjustment screw (**H**) as far as it will go clockwise (to the right).
  - Do not turn the screw beyond the natural right or left limit of rotation, as this will destroy the potentiometer.
  - Use the "Temp" turn dial (**E**) to adjust the set-point temperature to desired safety temperature ("Safe Temp").
  - Turn the "Safe Temp" set screw (**H**) slowly to the left until the heating function switches off and the signal light (**I** or **J**) goes off. Wait for about one more minute to ensure that the signal light (**F**) going off does not merely indicate the end of a switching cycle of the heater. If the signal light (**I** or **J**) turns on again within this time, you must repeat the procedure, slowly turning the "Safe Temp" (**H**) adjustment screw to the left again.
- When the signal light (**I** or **J**) does not light up again within one minute, the safety temperature is correctly set.
- After this, turn set "Safe Temp" adjustment screw (**H**) back slightly to the right (clockwise).
  - After you have turned the unit switch (**A**) of and back on again, the device is ready for operatin.

NOTE: When the "Safe Temp" (**H**) adjustment screw is set to the minimum level (as far as it will go to the left) the heating function is turned off.

Triggering of the safety circuit is shown in the following ways:  
On **RH basic 2** and **RH basic KT/C** versions: the warning light (I) will flash two times per second. (Note: regular flashing indicates that the safety circuit has been triggered; this warning light remains permanently lit to indicate that a standard heating cycle is in progress).  
On the **RH digital KT/C** version: the letter “-S-” will be shown on the LED display (G).

## Controlling the Medium Temperature via ETC 1 or Contact Thermometer

The control of the medium temperature via ETC 1 or contact-thermometer should be preferred. Its advantages are: just a short heating-up phase after setting the rated temperature, practically no temperature drift and only a minor temperature waviness.

The rear of the instrument houses the diode bushing for connection the ETS-D or contact thermometer, or for plugging the contact plug.

**The electronics of the instrument supply a test current that has to flow over plug pins 3 and 5 of the diode bushing in order for the heating plate to heat.**

### Safety contact thermometers

acc. to DIN 12 878 class 2 or acc. Gerstel are connected with a three-wire cable, the test current flows through the contact thermometer.

#### Safety function:

If the test current is interrupted because of e.g. breakage of contact thermometer or falling out of the cable plug, the heating cuts off.

Contact thermometer without safety circuit  
acc. to DIN 12 878 class 0. The instrument only heats if the test current circuit is closed by an electrical connection of the plug pins 3 and 5.

2-wire connecting cables:  
Connect plug pins 3 and 5 of the instrument plug.

3-wire connecting cables:  
Here the test current circuit can also be produced in the terminal head of the contact thermometer (connect plug pins 2 and 3).  
A 3-wire cable with the required bridge is available (accessories).

The **rated temperature** is set on the **contact thermometer** by means of the rotary magnetic head, whereby the upper edge of the traveling nut is the reference point on the rated value scale.

The required **surface temperature of the heating plat** is preselected on the **temp-rotary knob** of the instrument.

If the temp-rotary knob is set to the maximally adjustable temperature, the heating-up takes place as rapidly as possible, yet the medium temperature overshoots the rated value set on the contact thermometer.

If the temp-rotary knob is set to twice the rated value, the result is good compromise between rapid heating-up and overshooting. If the temp-rotary knob is set precisely to the rated temperature, the medium does not reach the rated temperature owing to a drop in temperature between heating plate and medium.

The **rated temperature** is set on the **ETS-D**.

If you are using an ETS-D4, we recommend using the ETS-D4 in On/Off mode (two-point control) at higher medium temperatures.

## Error - Messages

Version **RH digital KT/C:**

The LED displays (**C** and **D**) will start to flash and display a code when a fault is present. The fault codes and troubleshooting procedure are as follows:

<b>Fault code</b>	<b>LED</b>	<b>Cause</b>	<b>Solution</b>
<b>E 2</b>	G	Inner temperature too high	Switch off the unit and allow it cool down. Check that the ventilation slots on the bottom are not blocked and that cooling air can circulate freely.
<b>- S -</b>	G	Safety circuit is triggered	Reduce the temperature setpoint or reset the safety circuit where necessary.
<b>E 2</b>	D	No pulse to the reading fork (faulty reading fork, faulty motor or seized motor)	Switch off the unit. Trace the external cause blocking the stirring drive. If you cannot find and visible cause, clean the unit and return it for repair.
<b>E6</b>	G	Not put contact plug Break and/or short-circuit of the contact-thermometers	Contact plugs put  exchange the contact thermometer
Other codes			Switch off the unit. Clean and return for repair.

Version **RH basic 2** and **RH basic KT/C:**

The warning light (**I**) flashes to indicate that a fault is present. To remove a fault condition, proceed as follows:

- Switch off the unit and allow it to cool down.
- Trace and remove the cause of the fault.
- Switch on the unit.

If the fault persists, clean the unit and return it for repair.

## Mounting - Support rod H 16 V

The H 16V support rod can be screwed into the threading on the right behind the housing as follows:

- Remove the plastic cap and washer from the support rod and screw off the hex nut.
- Place the washer and hex nut on the beginning of the threading on the housing and hold them in place.
- Screw the support rod into the nut and threading of the housing and counter-tighten the hex nut.

## Maintenance and Cleaning

The IKA RH basic 2; RH basic KT/C or RH digital KT/C is maintenance-free. It is subject only to the natural wear and tear of components and their statistical failure rate.

**When ordering spare parts, please give the manufacturing number shown on the type plate, the machine type and the name of the spare part.**

Please send in equipment for repair only after it has been cleaned and is free from any materials which may constitute a health hazard. Use only cleansing agents which have been approved by IKA to clean IKA devices. To remove use:

Dyes	isopropyl alcohol
Construction materials	water containing tenside / isopropyl alcohol
Cosmetics	water containing tenside / isopropyl alcohol
Foodstuffs	water containing tenside
Fuels	water containing tenside

For materials which are not listed, please request information from IKA. Wear the proper protective gloves during cleaning of the devices. Electrical devices may not be placed in the cleansing agent for the purpose of cleaning.

Before using another than the recommended method for cleaning or decontamination, the user must ascertain with the manufacturer that this method does not destroy the instrument.

Version **RH basic KT/C** and **RH digital KT/C**

#### **Information for Care and Maintenance of the HeatingPlate with Technical Enamel Coating**

The technical enamel makes the heating plate easier to care for and more resistant to acids and bases. Because of it, however, the heating plate is also more susceptible to extreme fluctuations in temperature and the force of impact. This can result in cracks forming or the coating flaking off.

Make certain that the bottom of the placing vessel is even, clean and dry. The bottom of the placing vessel must not have any sharp grooves, sides or edges. Remove residues of bases and immediately. We recommend most strongly that you clean the heating plate regularly.

## Accessories

**RS 1** set of stirring bars

**H 15** bath top

Vers. KT

**H 36** holding rod

**H 44** cross sleeve

**H16V** support rod

**RSE** PTFE stirring bar remover

**ETC 1** IKATRON

**ETS-D** IKATRON

#### **Recommended IKA-stirres**

**TRIKA**-stirrer PTFE, triangular, length 55mm

**IKAFLO**N-stirrers PTFE

dia 8mm lengths 25, 30, 40 and 50 mm

## Technical data

		<b>RH basic 2</b>	<b>RH basic KT/C</b>	<b>RH digital KT/C</b>
design voltage:	<b>VAC</b>	220 - 240 ±10%	220 - 240 ±10%	220 - 240 ±10%
	or <b>VAC</b>	115 ±10%	115 ±10%	115 ±10%
	or <b>VAC</b>	100 ±10%	100 ±10%	100 ±10%
design frequency:	<b>Hz</b>	50/60	50/60	50/60
motor power consumption:	<b>W</b>	15	15	15
motor power output:	<b>W</b>	2	2	2
speed range:	<b>rpm</b>	100 to 2000	100 to 2000	100 to 2000
adjustment of the speed control:		infinitely variable setting	infinitely variable setting	infinitely variable setting
display of the speed control:		rotary knob on the front	rotary knob on the front	rotary knob on the front
motor:		scala 0 - 6	scala 0 - 6	LED-display four digit (0 - 2000 rpm)
max. stirring volume (water):	<b>ltr</b>	shaped pole motor	shaped pole motor	shaped pole motor
heating plate:		10	15	15
heating plate dimensions:	<b>mm</b>	1.4301	1.4301 with enamel coating	1.4301 with enamel coating
heating power:	<b>W</b>	Ø 125	130 x 130	130 x 130
	<b>W</b>	400 (230VAC)	500 (230VAC)	500 (230VAC)
	<b>W</b>	400 (115VAC)	500 (115VAC)	500 (115VAC)
	<b>W</b>	360 (100VAC)	450 (100VAC)	450 (100VAC)
temperature range:	<b>°C</b>	RT ..... 320	RT ..... 320	RT ..... 320
temperature setting:		rotary knob on the front	rotary knob on the front	rotary knob on the front
temperature display:		scala 0 - 6	scala 0 - 6	LED-display four digit (0°C - +320°C)
contact thermometer connection:		-----	DIN 12 878	DIN 12 878
infinitely variable safe temperature:	<b>°C</b>	-----	100 ....400	100 ....400
fixed adjustment safe temperature:	<b>°C</b>	400	-----	-----
perm. ambient temperature:	<b>°C</b>	+5 to +40	+5 to +40	+5 to +40
perm. humidity:	<b>%</b>	80	80	80
perm. on time:	<b>%</b>	100	100	100
screened for radio interference:		under VDE 0871	under VDE 0871	under VDE 0871
protection to DIN 60529:		IP 21	IP 21	IP 21
contamination level:		2	2	2
overvoltage categorie:		II	II	II
protection class :		1(protective earth)	1(protective earth)	1(protective earth)
fuse:	<b>A</b>	10AT Id-Nr. 2755400	10AT Id-Nr. 2755400	10AT Id-Nr. 2755400
operation at a terrestrial altitude:	<b>m</b>	max. 2000 above sea level	max. 2000 above sea level	max. 2000 above sea level
dimensions:	<b>mm</b>	168 x 220 x 105	168 x 220 x 105	168 x 220 x 105
weight:	<b>kg</b>	2,4	2,4	2,4

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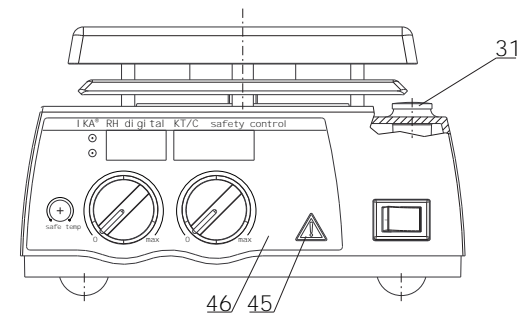
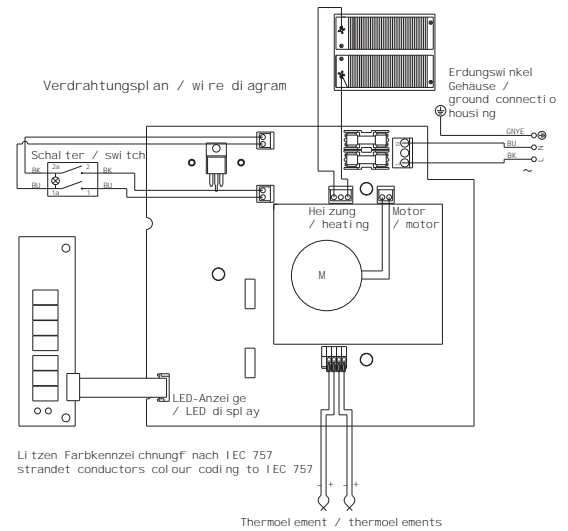
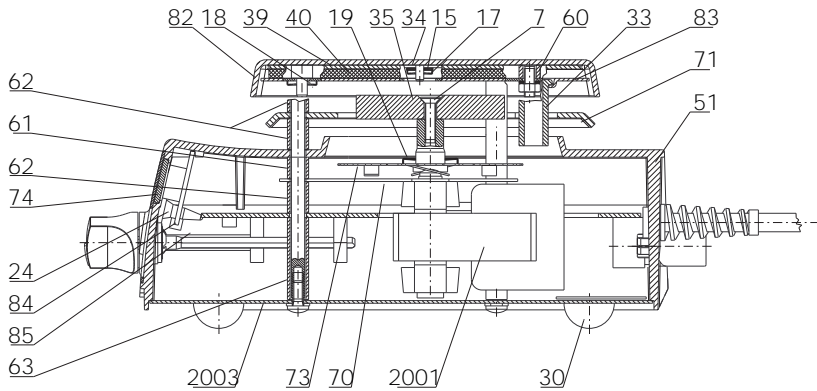
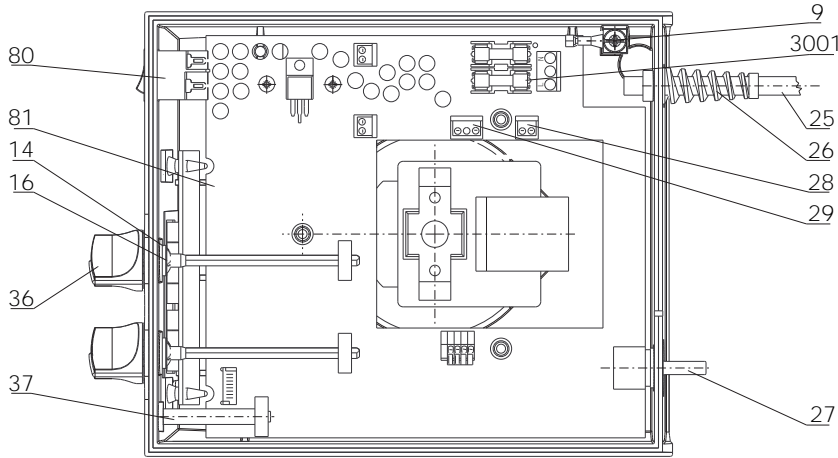
## List of spare parts

Item.	Designation
7	Countersunk screw M4 x16
9	Screw with washer assembly
14	Plastic washer
15	Washer A3,2 (Vers. KT/C)
16	Spring nut
17	Spring nut (Vers. KT/C)
18	Spring nut (Vers. KT/C)
19	Spring nut (Vers. KT/C)
24	Distance holder LP (Vers. digital KT)
25	Connection line
26	Mains lead cleat
27	Cap (Vers. RH basic 2)
27	Contact plug (Vers. KT/C)
28	Connector 2pol
29	Connector 3pol
30	Base
31	Screwed plug
33	Ceramic bushing (Vers. KT/C)
34	Thermoelement (Vers. KT/C)
35	Magnet
36	Rotary knob
37	Potentiometer arbor (Vers. KT/C)
39	Tubular heating element (Vers. RH basic 2)
39	Foil heating (Vers. KT/C)
40	Insulating (Vers. KT/C)
45	Front foil
46	Danger sign
49	Glass silk hose (Vers. RH basic 2)
51	Housing upper part
59	Bushing 85 Flat-strip lead compl. (Vers. digital)
60	Bushing 2001 Motor compl.
61	Bushing 2002 Thermoelement compl.
62	Bushing 2003 Bottom plate compl.
63	Bushing 3001 Fuse 6,3A
70	Support plate
71	Protection - sheet metal
73	Toothed washer
74	Window employment (Vers. basic 2)
74	Filter foil (Vers. digital)
75	Heat dissipator plate (Vers. RH basic 2)
76	Pressing sheet metal (Vers. RH basic 2)
77	Placing plate (Vers. RH basic 2)
80	Rocker switch
81	PCB control
82	Placing plate (Vers. KT/C)
83	Pressing sheet metal (Vers. KT/C)
84	PCB display (Vers. digital)

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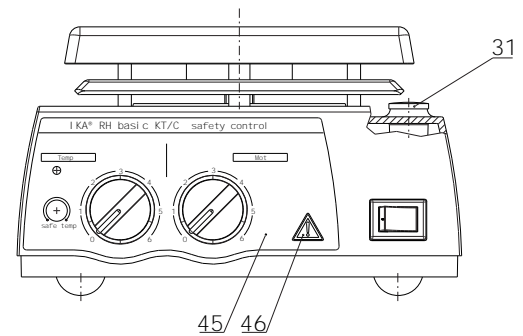
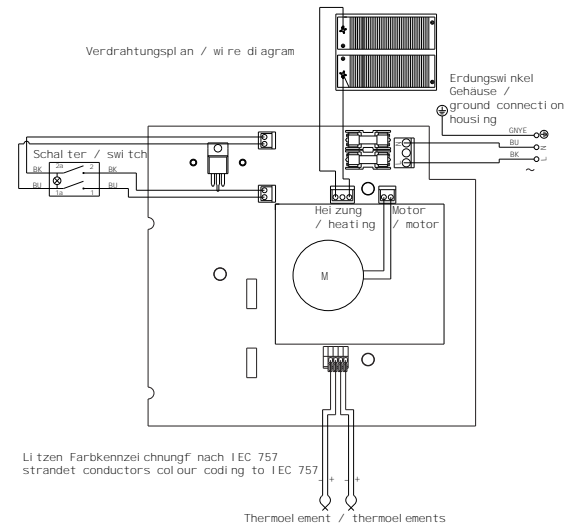
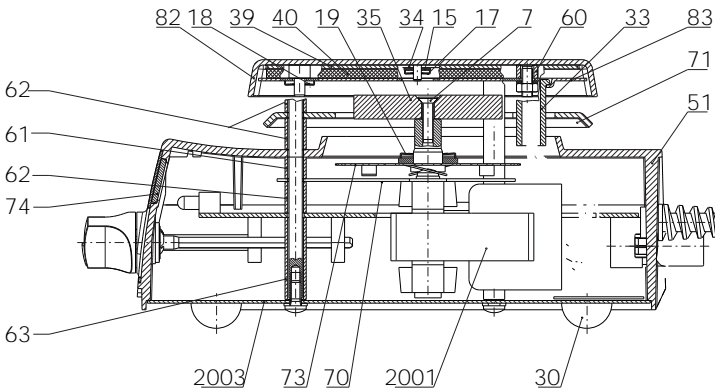
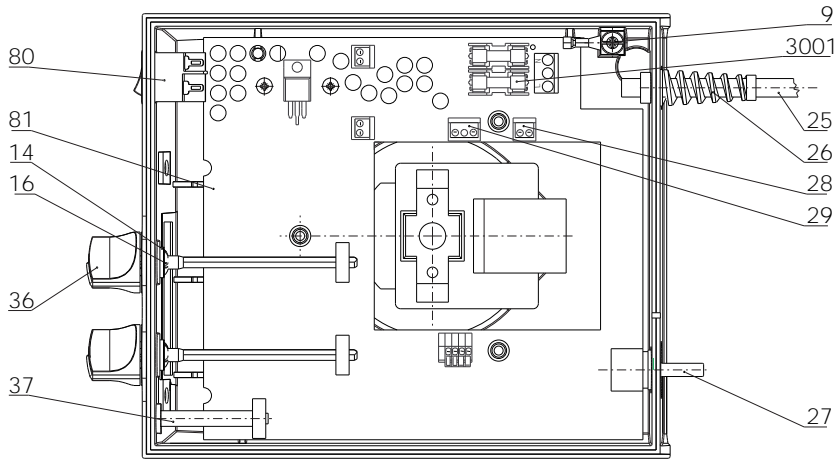
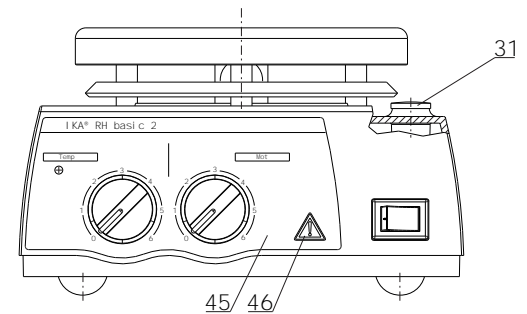
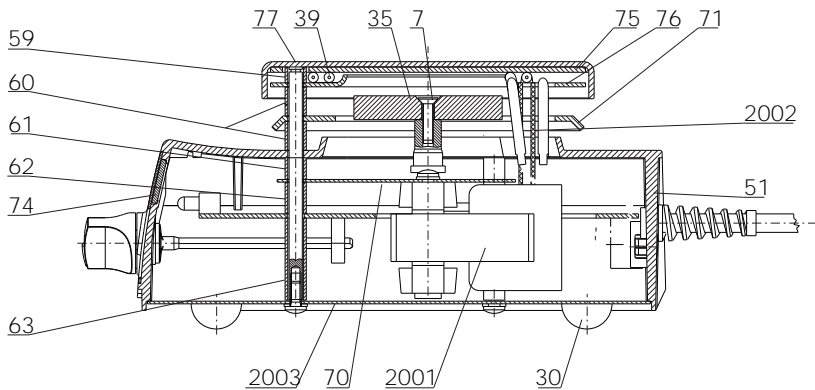
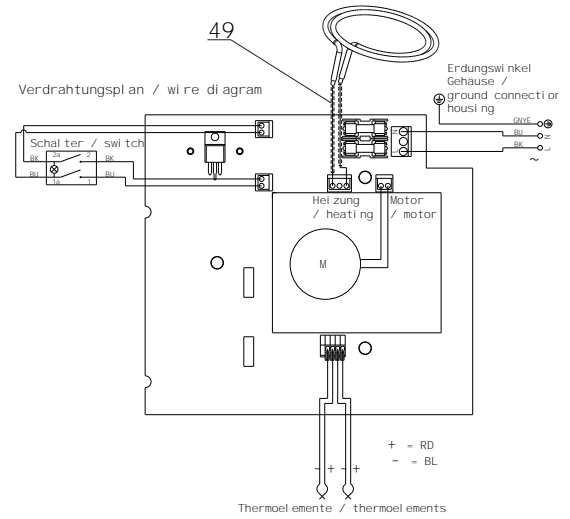
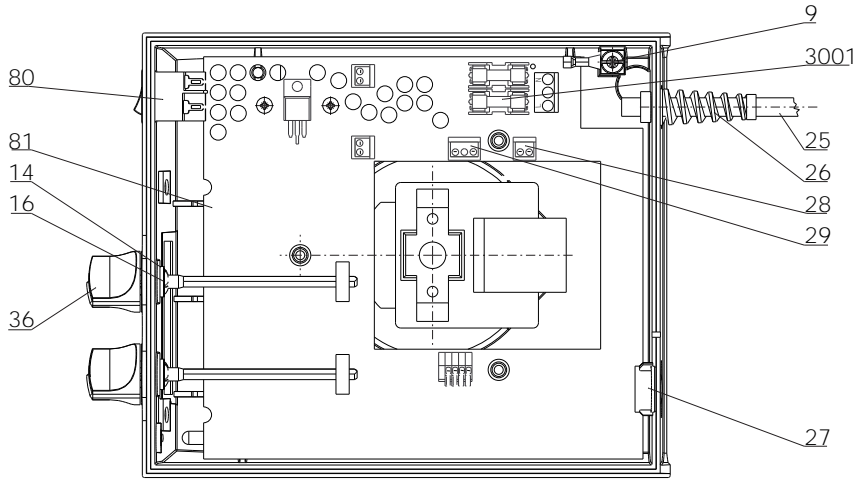
China

---

IKA® Works Guangzhou  
173 - 175 Friendship Road  
Guangzhou  
Economic and Technological  
Development District  
510730 Guangzhou, China  
Tel. +86 20 8222-6771  
Fax +86 20 8222-6776  
sales@ikagz.com.cn

Japan

---

IKA® Japan K.K.  
3-5-8 Yokonuma-cho,  
Higashiosaka-city, Osaka  
577-0808 Japan  
Tel. +81-6-6730-6781  
Fax +81-6-6730-6782  
info@ika.ne.jp

India

---

IKA® India Private Limited  
814/475, Survey No. 129/1  
Mysore Road, Kengeri  
Bangalore - 560 060  
Karnataka, India  
Tel. +91 80 26253900  
Fax +91 80 26253901  
info@ika.in

Asia  
Australia

---

IKA® Works (Asia)  
Sdn Bhd  
No. 17 & 19, Jalan PJU 3/50  
Sunway Damansara  
Technology Park  
47810 Petaling Jaya  
Selangor, Malaysia  
Tel. +60 3 7804-3322  
Fax +60 3 7804-8940  
sales@ika.com.my

Korea

---

IKA® Korea Co LTD  
1710 Anyang Trade Center  
1107 Buhung-dong,  
Dongan-gu  
Anyang City, Kyeonggi-do  
Post code: 431-817  
South Korea  
Tel. +82 31-380-6877  
Fax +82 31-380-6878  
michael@ikakorea.co.kr

Brasilia

---

IKA® Works Inc.  
Rua São Bento, 701 Sala 1  
CEP 13160-000 Centro -  
Artur Nogueira - SP  
Brasil, South America  
Tel. +55 19 3877 2399  
Fax +55 19 3877 2399  
fcabra@ika.net