IKAMAG® KMO 2 basic

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OPERATING INSTRUCTIONS EN 7

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DE

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CE-DECLARATION OF CONFORMITY

EN

We declare under our sole responsibility that this product corresponds to the regulations 89/336EEC and 73/023EEC and conforms with the standards or standardized documents EN 60 010-1; EN 50 081; EN 50 082; EN 55 014 and EN 60 555.

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Guarantee

You have purchased an original IKA laboratory machine which meets the highest engineering and quality standards.

In accordance with IKA guarantee conditions, the guarantee period is 24 months. For claims under the guarantee please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

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Contents

	Page
Guarantee	2
Safety instructions	7
Correct use	7
Unpacking	7
Commissioning	8
General information	8
Function Stirring	8
General	8
Maintenance and cleaning	8
Associated standards and regulations	9
Accessories	9
Technical data	9
List of spare parts KMO 2 basic	14
Spare parts diagram KMO 2 basic	15

Safety instructions

The instrument may not be operated in rooms with explosion hazards.

Furthermore, it has to be placed upon a fire-proof and/or not burnable surface.

Attention: Magnetism! Effects of the magnetic field have to be taken into account (e.g. data carriers, cardiac, cardiac pacemakers...).

When using PTFE-coated magnetic bars, please pay attention to the following: Chemical reactions of PTFE occur in contact with molten or dissolved alkaline and alkaline earth metals, as well as with fine-particled powders of metals of the 2. and 3. group of the periodical system at temperatures above 300 - 400°C. Only elementary fluorine, chlorine trifluoride and alkaline metals do attract PTFE, halogen hydrocarbons have a reversibly swelling effect.

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Risk of burns!

The motor mounting screws on the underside of the device can reach temperatures in excess of 100°C! The shape of the housing at the side protects the screws against accidental contact.

When carrying or lifting the device, take care not to touch the heads of the screws with your fingers!

Even in the case of repair work, the unit must only be opened by a qualified engineer. The mains plug must be pulled before opening the unit. The voltage-carrying components inside the unit may continue to carry a voltage for some time after the mains plug is pulled.

Correct use

The IKAMAG KMO 2 basic is a small, unheated magnetic stirrer. The device is suitable for mixing up to 3 liters of low-viscosity media in receptacles placed on the device. The built-in stirrer drive allows the substances to be stirred by means of a magnetic rod in the receptacle (40 and 50 mm recommended). The intensity of mixing is dependent on the speed of the motor (approx. 100 to 1100 rpm) and the size of the magnetic rod.

Unpacking

Please unpack the equipment carefully and check for any damages. It is important that any damages which may have arisen during transport are ascertained when unpacking. If applicable a fact report must be set immediately (post, rail or forwarder).

The delivery scope covers: an IKAMAG KMO 2 basic and operating instructions.

Commissioning

Check whether the voltage specified on the type plate matches the mains voltage available. The power socket used must be earthed (protective earth conductor contact). If these conditions are met, the device is ready to operate after plugging in the mains plug. If these procedures are not followed, safe operation cannot be guaranteed and/or the equipment may be damaged.

Observe the ambient conditions (temperature, humidity, etc.) listed under Technical Data.

General information

A split-pole motor carries a powerful permanent magnet and a shielding plate. This shielding plate redirects the magnetic field lines upwards, thus amplifying the magnetic field in the direction of the baseplate. It also carries the ring gear used to measure the speed of rotation. The optoelectronic controller receives a pulse from each tooth on the shielding plate. The time interval between each pulse is a measure of the actual rotational speed. A downstream comparator calculates whether the selected speed setting matches the actual speed and opens the power supply to the motor whenever the actual speed drops below the setting.

Function Stirring

The controls are located on the front panel. Press the power switch to activate the device. When switched on the green indicator lamp in the power switch is lit. After switching on the machine starts to run at the most recent speed setting. The speed of the motor can be set on the rotary knob with scale.

General

An M10 threaded bushing is located on the left at the rear of the baseplate. This can be used to attach a stand rod for fastening measuring and metering equipment.

CAUTION! The device must not be lifted, carried or fastened to a stand by the stand rod in the threaded bushing. The device is intended exclusively for use on a hard, level work surface.

NOTE! The device will become warm during operation due to power losses from the motor and the electronics. In order to dissipate this heat and to avoid heating the baseplate to a critical degree, fan blades are integrated into the motor to circulate the air inside the device. For this reason, never obstruct the ventilation slots at the bottom and rear of the device, and ensure that an adequate distance is left between the device and other objects.

Maintenance and cleaning

The stirrer RW 11 basic are maintenance-free. It is subject only to the natural wear and tear of components and their statistical failure rate.

When ordering spare parts, please give the manufacturing number shown on the type plate, the machine type and the name of the spare part.

Please send in equipment for repair only after it has been cleaned and is free from any materials which may constitute a health hazard. Use only cleansing agents which have been approved by IKA to clean IKA devices.

To remove	use:
Dyes	isopropyl alcohol
Construction materials	water containing tenside / isopropyl alcohol
Cosmetics	water containing tenside / isopropyl alcohol
Foodstuffs	water containing tenside
Fuels	water containing tenside

For materials which are not listed, please request information from IKA.
Wear the proper protective gloves during cleaning of the devices.
Electrical devices may not be placed in the cleansing agent for the purpose of cleaning.

Associated standards and regulations

Construction in accordance with the following safety standards
EN 61 010-1 UL 3101-1 CAN/CSA C22.2 (1010-1)

Construction in accordance with the following EMC standards
EN 50 082-1 EN 55 014 EN 60 555-2; -3

Associated EU guidelines
EMC-guidelines: 89/336/EWG
Machine guidelines: 73/023/EWG

Accessories

RS 1	Set of stirrers	H16V	Support rod
RSE	Stirrer remover	H 15	Bath attachment
DZM control.m	Rpm counter		

Recommended IKA stirrers

TRIKA stirrer, PTFE, triangular, 42 and 55 mm long

IKAFILON stirrers, PTFE, 7 mm dia.

lengths 30, 40, 50 and 60 mm

Technical data

Design voltage:	VAC	230±10%	
	or	VAC	115±10%
Design frequency:	Hz	50/60	
Power consumption:	W	14	
Power output:	W	4	
Speed range:	rpm	ca. 100 to 1100 infinitely variable	
Max. stirring quantity (Water)	ltr	3	
Perm. ambient temperature:	°C	+5 to +40	
Perm. relative humidity:	%	80	
Perm. ON duration:	%	100	
Radio-screened:		acc. to VDE 0871	
Protection to DIN 40 050:		IP 21	
Housing:		PA 6.6 M30	
Baseplate:		1.4301	
Dimensions:	mm	140 x 200 x 75	
Weight:	kg	1,4	

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List of spare parts

KMO 2 basic

Item	Designation
9	connection line
10	strain relief
11	protective cap
14	pedestal
21	control knob
22	potentiometer shaft
23	clamping bushing
26	front panel
27	Zierkappe
30	housing - upper part
31	housing - lower part
35	printed circuit board
37	hood
40	rocker switch
2001	drive unit, assembly

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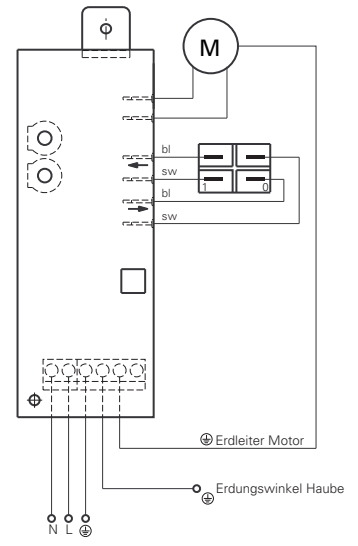
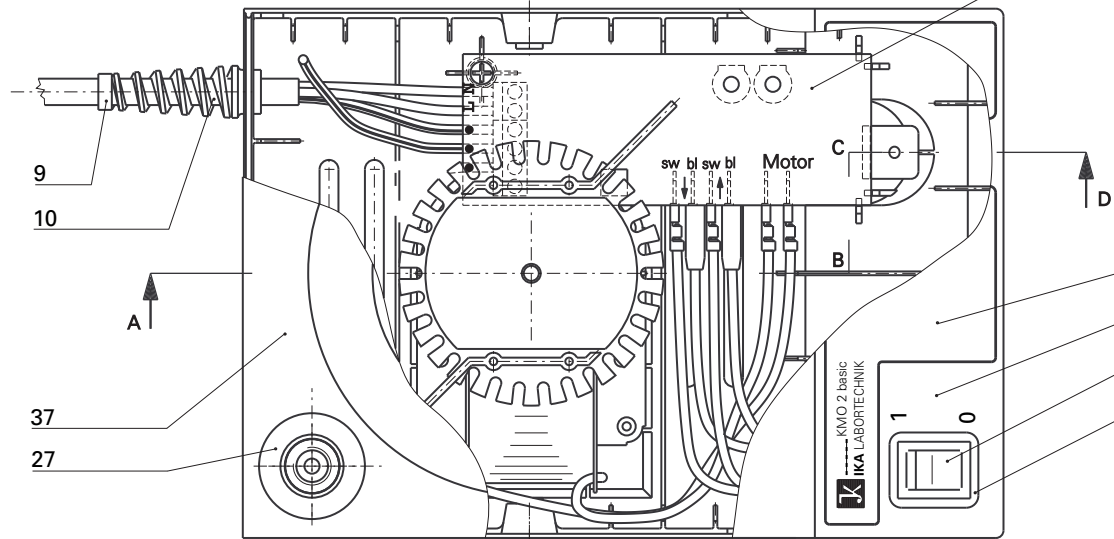
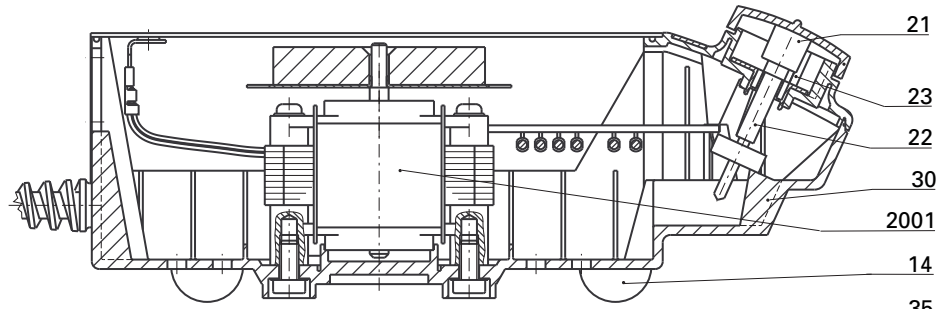
India

IKA® - Werke
GmbH & Co.KG
Liaison Office India
No. 31 (Old No. 264)
1st Floor, 10th Cross
1st "N" Block, Rajajinagar
560 010 Bangalore
Tel.: +91 80-41157736
Fax: +91 80-41157735
E-Mail: info@ikaindia.com

Asia
Australia

IKA® Works (Asia)
Sdn Bhd
No. 17 & 19, Jalan PJU 3/50
Sunway Damansara
Technology Park
47810 Petaling Jaya
Selangor, Malaysia
Tel.: +60 3 7804-3322
Fax: +60 3 7804-8940
E-Mail: sales@ika.com.my

Korea

IKA® Korea Co LTD
1710 Anyang Trade Center
1107 Buhung-dong,
Dongan-gu
Anyang City, Kyeonggi-do
Post code: 431-817
South Korea
Tel.: +82 31-380-6877
Fax: +82 31-380-6878
E-Mail: michael@ikakorea.co.kr

Brasilia

IKA® Works Inc.
Av. das Américas,
15700, sala 235
Recreio dos Bandeirantes
CEP 22790-701
Rio de Janeiro, RJ
Brasil
Tel.: +55 21 2487-7743
Fax: +55 21 2487-7743
E-Mail: fcabral@ika.net